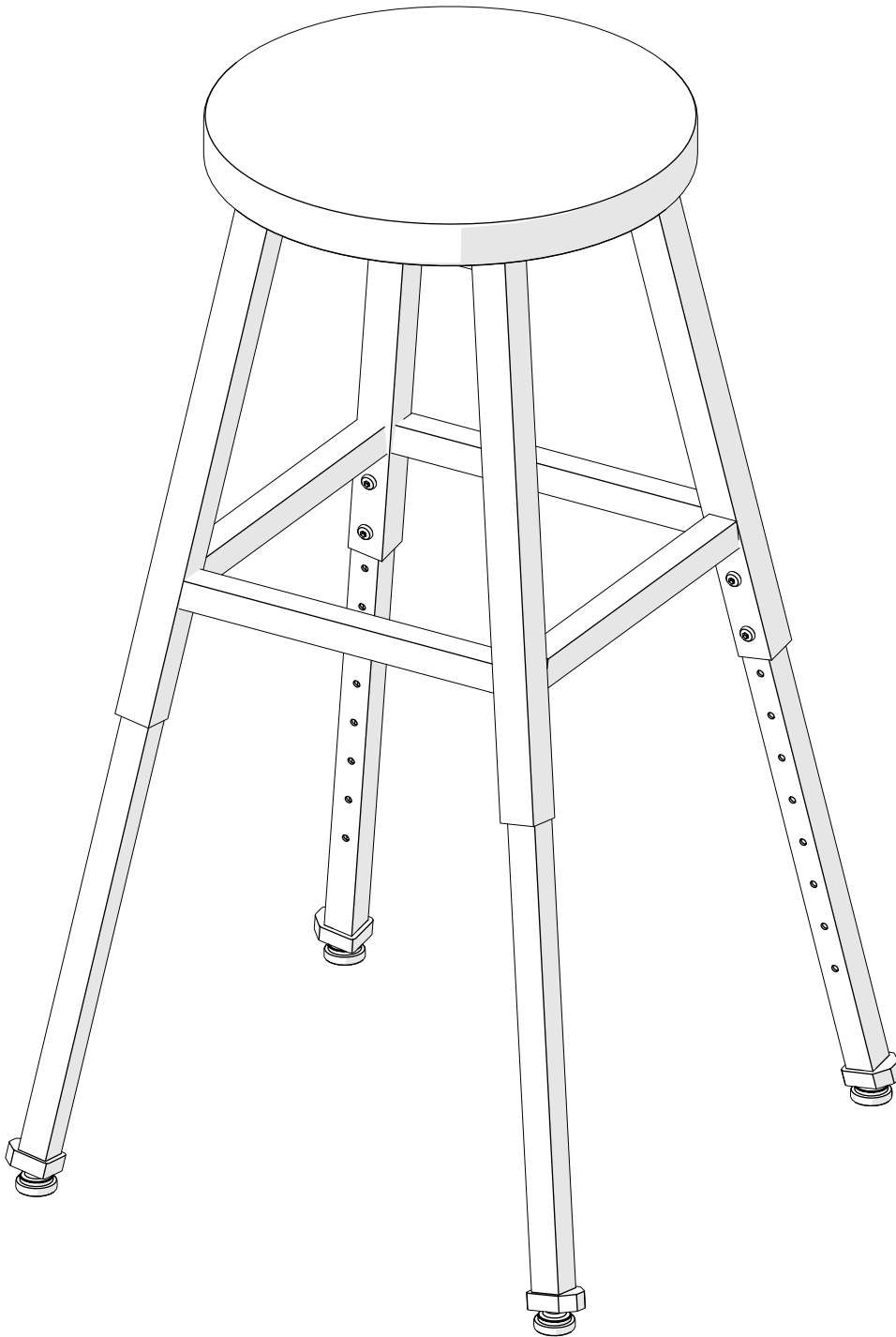
# Lab Stool

34419R

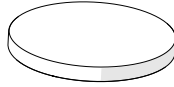
 **ASSEMBLY INSTRUCTIONS**  **DO NOT THROW AWAY!**

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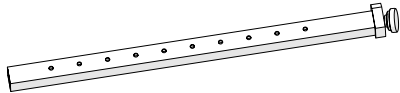
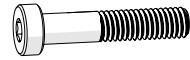
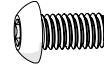
x8

## WARNINGS AND DISCLAIMERS

Review full assembly instructions and check that all parts are present before beginning assembly. Heavy or large furniture may require multiple persons in order to properly assemble. If you have questions regarding assembly, contact the manufacturer before continuing. Damage due to incorrect assembly may not be covered by warranty. Do not use product unless all screws are tight. Check all screws for tightness at least every six months. If parts are broken, missing, damaged, or worn, stop use of the product until repairs are made by your dealer using factory authorized parts.

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## WARRANTY

For full product warranty details, please visit:

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<http://moorecoinc.com/warranty>

## CARE & MAINTENANCE

For questions on assembly or missing/damaged parts, please contact us using the following information.

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Central Time)

## PRODUCT REGISTRATION

To register your product for warranty, please visit:

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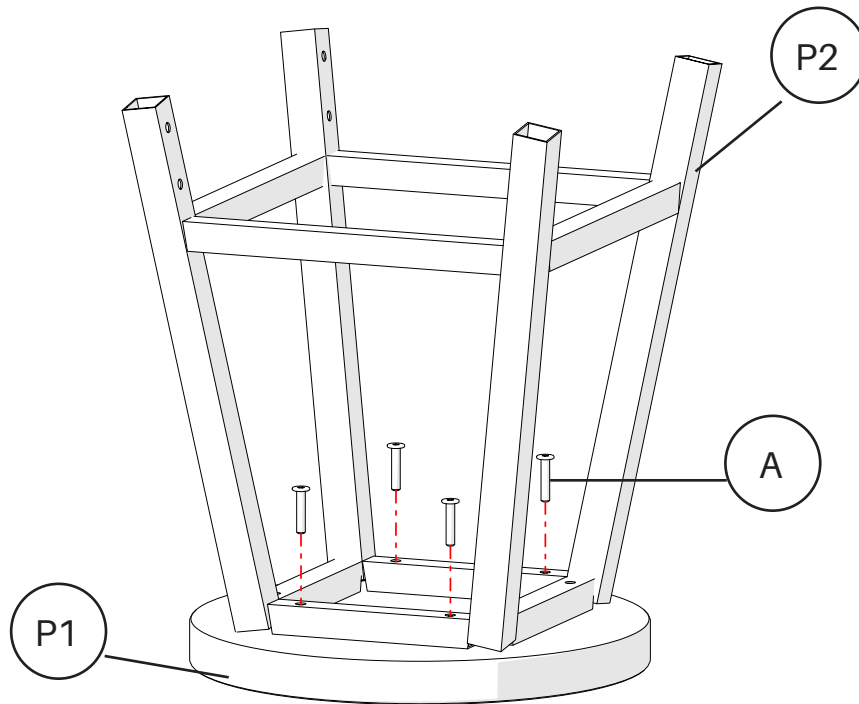
<https://moorecoinc.com/register>


## STEP 1

Place the seat (P1) upside-down on a non-scratch surface. Attach the frame (P2) to the seat by screwing the M6 x 30mm (A) screws through the holes in the frame into the inserts on the seat.

## STEP 2

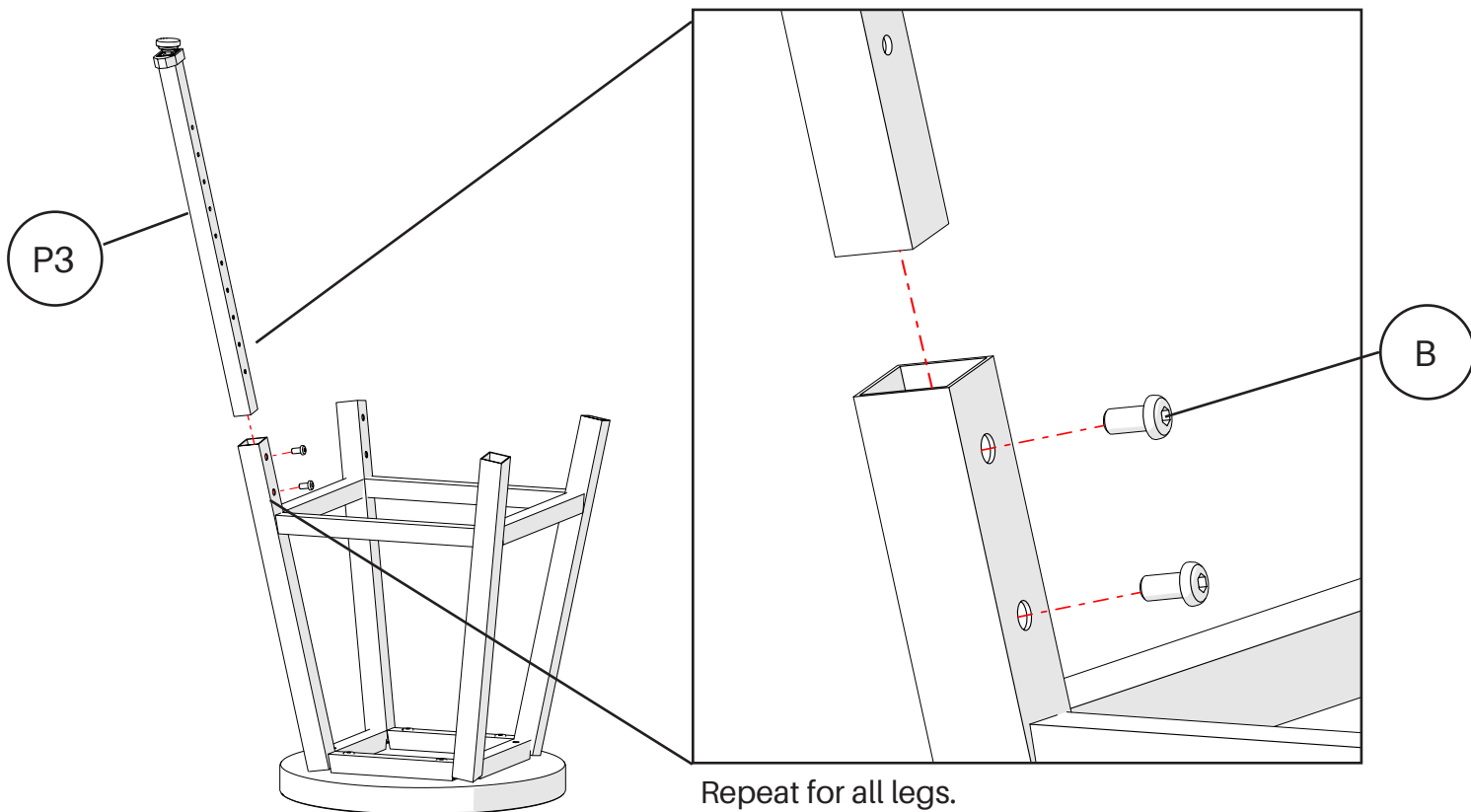
Attach the adjustable leg (P3) to the frame using the M6 x 12 mm (B) button head screws. Repeat for all legs.

---

**STEP 4**

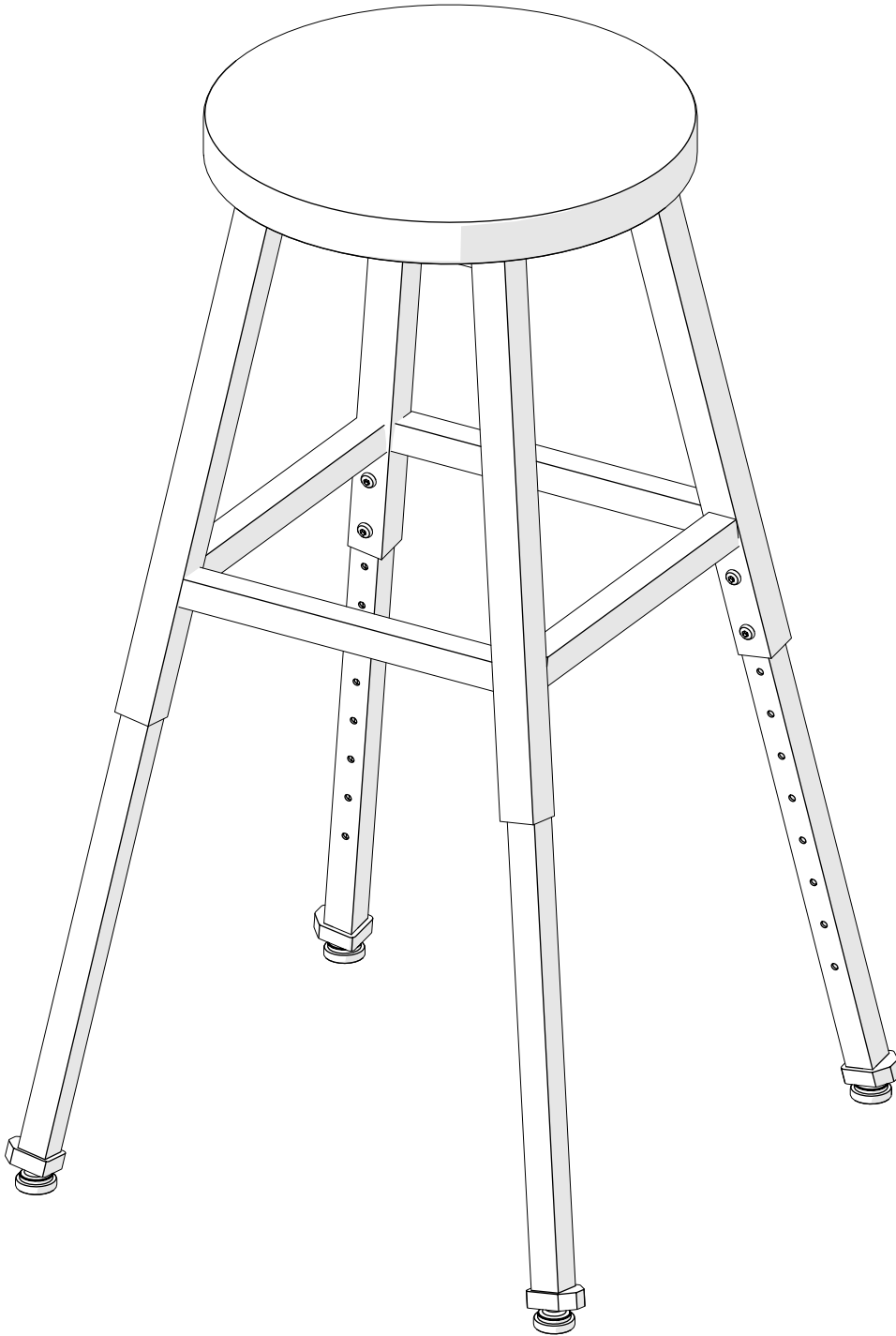Flip stool over onto its feet.

---

**THIS COMPLETES ASSEMBLY INSTRUCTION**